

TeSR™2, a Xeno-Free Version of mTeSR™1 for Maintenance of Human ES and iPS Cells

TeSR™2 is a xeno-free medium for long-term maintenance of human embryonic stem (ES) and induced pluripotent stem (iPS) cells. The TeSR™ family of feeder-free maintenance media includes mTeSR™1, TeSR™2 and TeSR™-E8™, which are based on published formulations from the laboratory of James Thomson.¹⁻³ Closely related to mTeSR™1, the most widely published feeder-free human pluripotent stem cell (hPSC) medium, TeSR™2 has a modified formulation, but with all xeno-free components, to produce a more defined medium.²

As a xeno-free alternative, TeSR™2 provides the same high-quality and robust system for basic research, stem cell banking, scale-up studies and pre-clinical applications. TeSR™2, can also be used with Vitronectin XF™ (Catalog #07180) as a substrate to culture cells in a completely xeno-free system.

Comparable with mTeSR™1

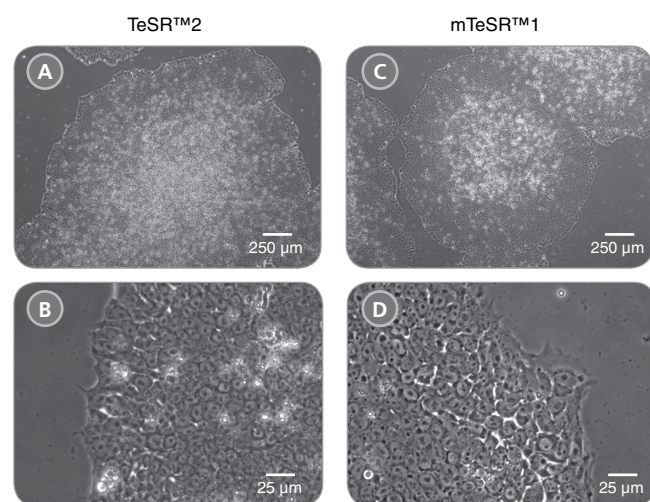


FIGURE 1. Morphology of hPSCs Maintained in TeSR™2 is Comparable to hPSCs Cultured in mTeSR™1

(A,B) Undifferentiated human ES (H9) cells cultured on Corning® Matrigel® matrix in TeSR™2 retain the prominent nucleoli and high nuclear-to-cytoplasmic ratio characteristic of this cell type. Densely packed cells and multilayering are apparent when cells are ready to be passaged. **(C,D)** H9 cells cultured under the same conditions in mTeSR™1 exhibit comparable morphology.

Advantages:

XENO-FREE. TeSR™2 is a more defined version of mTeSR™1, free of xenogenic components.

COMPATIBLE. Use with published mTeSR™1 protocols for a wide variety of applications.

ROBUST. Formulation contains recombinant human albumin to aid in lipid/nutrient transport and to protect cultures from cellular toxins and stresses.

INTEGRATED WORKFLOW. Maintain newly generated human iPS cells (reprogrammed using TeSR™-E7™ or ReproTeSR™) and hPSCs prior to directed downstream differentiation with STEMdiff™ products.

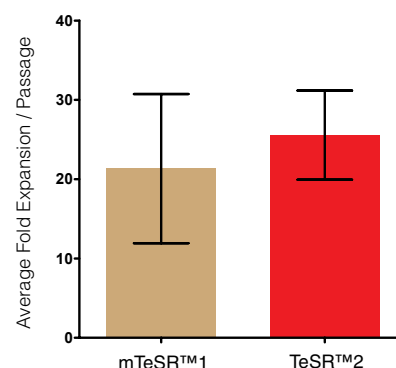


FIGURE 2. Fold and Cumulative Aggregate Expansion in TeSR™2

Graph shows the average fold expansion per passage \pm SEM obtained for human ES and iPS cells cultured in mTeSR™1 (brown) or TeSR™2 (red) with Corning® Matrigel® over 10 passages. Expansion was determined by counting the cell aggregates obtained at harvest and dividing by the number of cell aggregates seeded.

Note: This data is representative of cultures passaged after 5-6 days in culture; lower expansion should be expected if using shorter culture times.

TeSR™2

Defined, Feeder-Free and Xeno-Free hES and hiPS Cell Medium

Standardized Feeder-Free Maintenance

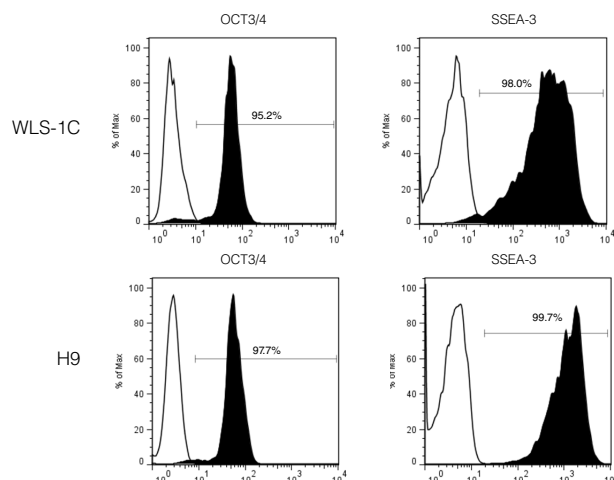


FIGURE 3. Human Pluripotent Stem Cells Cultured in TeSR™2 Retain Expression of Undifferentiated Cell Markers

Histogram analysis for H9 human ES and WLS-1C human iPS cells characterized using flow cytometry for undifferentiated cell markers (SSEA-3 and OCT3/4) after passaging in TeSR™2 for 21 passages (WLS-1C) and 18 passages (H9), respectively (filled histogram = sample, hollow histogram = secondary antibody only).



FIGURE 4. Human ES Cells Cultured Long-Term in TeSR™2 Retain Normal Karyotype

Chromosomal analysis of H9 hES cells cultured in TeSR™2 for 12 passages shows that normal karyotype is retained during passaging.

Product Information

PRODUCT	SIZE	CATALOG #
TeSR™2	1 Kit	05860
	10 Kits	05880
Vitronectin XF™	2 mL	07180
	1 Kit	07190

For a complete list of related products, including specialized cell culture and storage media, matrices, antibodies, cytokines and small molecules, visit www.stemcell.com/hPSCworkflow or contact us at techsupport@stemcell.com. Recommended antibodies include Anti-Human OCT4 (OCT3) Antibody, Clone 3A2A20 (Catalog #60093) and Anti-Mouse SSEA-3 Antibody, Clone MC-631 (Catalog #60061).

Copyright © 2016 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, Scientists Helping Scientists STEMdiff and ReLeSR are trademarks of STEMCELL Technologies Canada Inc. Corning and Matrigel are registered trademarks of Corning Inc. mTeSR, TeSR, E8, and E7 are trademarks of WARF. All other trademarks are property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.

STEMCELL TECHNOLOGIES INC.'S QUALITY MANAGEMENT SYSTEM IS CERTIFIED TO ISO 13485. PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED.

TeSR™2 Integrates Upstream or Downstream of Your Workflow

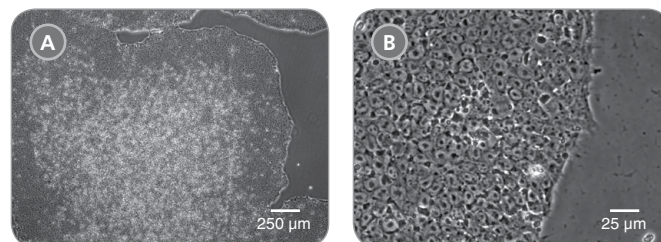


FIGURE 5. iPS Cell Colonies Generated with ReproTeSR™ Can Be Expanded in TeSR™2

(A,B) Representative images of iPS cell colonies generated using ReproTeSR™ and cultured in TeSR™2. Generated iPS cells retain the prominent nucleoli and high nuclear-to-cytoplasmic ratio characteristic of this cell type. Densely packed cells and multilayering are apparent when cells are ready to be passaged.

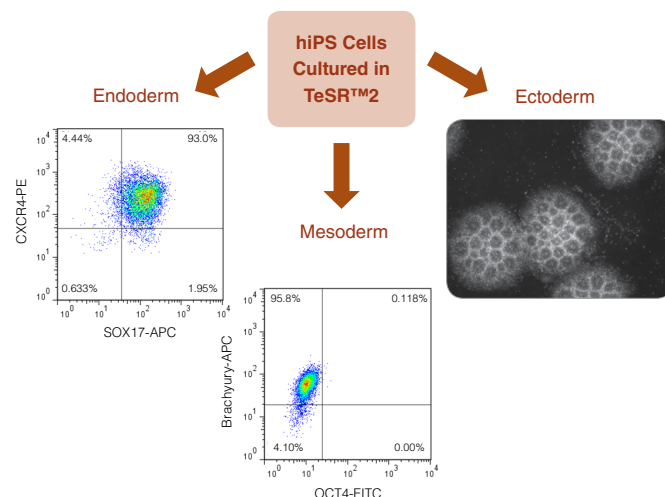


FIGURE 6. Directed Differentiation of TeSR™2-Maintained hiPS Cells

WLS-1C human iPS cells maintained in TeSR™2 were differentiated into all three germ layers. Endoderm specification was achieved using the STEMdiff™ Definitive Endoderm Kit (Catalog #05110). Mesoderm specification was demonstrated using a protocol modified from Lian X, et al.⁴ Ectoderm specification was demonstrated using STEMdiff™ Neural Induction Medium (Catalog #05835) to generate neural rosettes, a morphological hallmark of neural induction.

References

- Ludwig TE et al. (2006) Nat Methods 3(8): 637–46.
- Ludwig TE et al. (2006) Nat Biotechnol 24(2): 185–7.
- Chen G et al. (2011) Nat Methods 8(5): 424–9.
- Lian X et al. (2012) PNAS 109(27): E1848–57.