

Agar Leukocyte Conditioned Medium



Scientists Helping Scientists™ | WWW.STEMCELL.COM

TOLL FREE PHONE 1 800 667 0322 • PHONE +1 604 877 0713

INFO@STEMCELL.COM • TECHSUPPORT@STEMCELL.COM

FOR GLOBAL CONTACT DETAILS VISIT OUR WEBSITE

Source of colony-stimulating factors for assays of human hematopoietic progenitor cells

Catalog #02300

25 mL

Product Description

Agar Leukocyte Conditioned Medium (Agar-LCM) is a source of colony-stimulating factors for assays of human hematopoietic progenitor cells from light-density fractions of bone marrow and peripheral blood. It is suitable for the growth of colony-forming unit-erythroid (CFU-E), burst-forming unit-erythroid (BFU-E), colony-forming unit-macrophage (CFU-M), colony-forming unit-granulocyte (CFU-G), colony-forming unit-granulocyte, macrophage (CFU-GM), and colony-forming unit-granulocyte, erythroid, macrophage, megakaryocyte (CFU-GEMM) when added to MethoCult™ medium at a final concentration of 10% with the addition of erythropoietin (EPO; Catalog #78007).

Agar-LCM is prepared using normal human peripheral blood leukocytes in an agar-containing medium.

Properties

Storage: Store at -20°C.

Shelf Life: Stable until expiry date (EXP) on label.

Contains:

- Agar Leukocyte Conditioned Medium
- Fetal bovine serum (FBS; 10%)
- Iscove's Modified Dulbecco's Medium (IMDM)
- 1×10^{-4} 2-Mercaptoethanol (ME)

This product contains material derived from human plasma. Donors have been tested and found negative for HIV-1 and -2, hepatitis B, and hepatitis C prior to donation. However, this product should be considered potentially infectious and treated in accordance with universal handling precautions.

Directions for Use

Thaw Agar-LCM at room temperature (15 - 25°C) or overnight at 2 - 8°C.

NOTE: After thawing, some particulate matter may appear; the product may be clarified by centrifugation. There should be no loss in bioactivity as a result of this procedure. Keep on ice while in use.

PREPARATION OF METHOCULT™ H4230 (CATALOG #04230) CONTAINING AGAR-LCM FOR DETECTION OF HUMAN HEMATOPOIETIC PROGENITOR CELLS (BFU-E, CFU-GM, CFU-GEMM)

1. Thaw 80 mL bottle of MethoCult™ H4230 at room temperature or overnight at 2 - 8°C. Mix contents by vigorous shaking.
2. Add 10 mL of Agar-LCM. Add EPO to yield a final concentration of 3 U/mL. Add IMDM (e.g. Catalog #36150) to a total volume of 100 mL.
3. Mix medium by vigorous shaking for 1 - 2 minutes. Let stand for 5 minutes to allow bubbles to rise to the top.
4. Dispense 3 mL of medium into sterile culture tubes using a 6 or 12 mL syringe attached to a 16 Gauge Blunt-End Needle (Catalog #28110).

NOTE: If not used immediately, cap tubes tightly and store at -20°C. Thaw tubes at room temperature.

5. Prepare a suspension of human cells in IMDM with 2% FBS (e.g. Catalog #07700) at 10X the final concentration required for plating.
6. Add 0.3 mL of cells to 3 mL of MethoCult™ medium.

NOTE: This 1:10 (v/v) ratio of cells to medium gives the correct viscosity to ensure optimal CFU growth and morphology.

7. Vortex tubes vigorously for at least 4 seconds to mix contents thoroughly. Let stand for at least 5 minutes to allow bubbles to rise to the top.
8. Dispense 1.1 mL of MethoCult™ mixture containing cells into each of two 35 mm Culture Dishes (e.g. Catalog #27100) using a 3 mL syringe (e.g. Catalog #28230) attached to a 16 Gauge Blunt-End Needle.
9. Incubate for 12 - 14 days at 37°C in 5% CO₂ with ≥ 95% humidity.
10. Score colonies using an inverted microscope.

Notes and Tips

For additional information, refer to Technical Manual: Human Colony-Forming Unit (CFU) Assays Using MethoCult™ (Document # 10000005589), available at www.stemcell.com, or contact us to request a copy.

Related Products

For related products, including specialized culture and storage media, supplements, antibodies, cytokines, and small molecules, visit www.stemcell.com/HSPCworkflow. For available fresh and cryopreserved peripheral blood, cord blood, and bone marrow products, visit www.stemcell.com/primarycells.

PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED. FOR ADDITIONAL INFORMATION ON QUALITY AT STEMCELL, REFER TO WWW.STEMCELL.COM/COMPLIANCE.

Copyright © 2025 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, Scientists Helping Scientists, and MethoCult are trademarks of STEMCELL Technologies Canada Inc. All other trademarks are the property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.