

# Primary Cells

## Human Peripheral Blood Leukapheresis Pack, Fresh



Scientists Helping Scientists™ | [WWW.STEMCELL.COM](http://WWW.STEMCELL.COM)

TOLL FREE PHONE 1 800 667 0322 • PHONE +1 604 877 0713

[INFO@STEMCELL.COM](mailto:INFO@STEMCELL.COM) • [TECHSUPPORT@STEMCELL.COM](mailto:TECHSUPPORT@STEMCELL.COM)

FOR GLOBAL CONTACT DETAILS VISIT OUR WEBSITE

Catalog #	200-0092
	70500.2
	70500.1
	70500

Tenth Size
Quarter Size
Half Size
Full Size

## Product Description

Leukapheresis is performed on normal donors using Institutional Review Board (IRB)-approved consent forms and protocols. Approximately two to three blood volumes are processed using the Spectra Optia® Apheresis System (unless otherwise requested) to produce a full-sized Leukapheresis Pack.

Donor Status:	Normal
Characterization Criteria:	Cell count, viability, donor virus testing, age, sex, ethnicity, weight, height, smoking status, other information
Format:	Product is drawn directly into a sample collection bag containing anticoagulant.
Anticoagulant:	Acid citrate dextrose solution A (ACDA)

For donor details, refer to the lot-specific Certificate of Analysis.

## Stability and Storage

Product is shipped at ambient temperature and should be used or processed immediately upon receipt.

## Precautions

Donor Screening: Donors are screened for HIV-1, HIV-2, hepatitis B, and hepatitis C.

If the donor has been screened within 90 days of donation the product will be shipped with negative test results from donor screening.

If the donor has not been screened within 90 days of collection, a test sample will be taken at the time of donation and the product will be shipped before the screening results are available. In the unlikely event that a test result is positive, the customer will be contacted as soon as possible (usually within 24 - 72 hours from the time of shipment).

Donors have been tested and found to be negative for HIV-1, HIV-2, hepatitis B, and hepatitis C prior to donation. As testing cannot completely guarantee that the donor was virus-free, THIS PRODUCT SHOULD BE TREATED AS POTENTIALLY INFECTIOUS and only used following appropriate handling precautions such as those described in biological safety level 2. When handling this product, do not use sharps such as needles and syringes.

STEMCELL cannot guarantee the biological function or any other properties associated with performance of cells in a researcher's individual assay or culture systems. STEMCELL assures the cells will meet the specifications only when assessed, before washing, by our test methods.

FOR IN VITRO RESEARCH USE ONLY. NOT APPROVED FOR DIAGNOSTIC, THERAPEUTIC, OR CLINICAL APPLICATIONS.  
NOT APPROVED FOR HUMAN OR VETERINARY USE IN VIVO.

## Handling / Directions for Use

IMPORTANT: To determine the number of cells provided, a cell count must be done upon receipt and before any processing steps (e.g. washing). Cell loss is expected during wash steps and may be up to 30%. Use sterile technique when processing cells.

Remove a 20 µL aliquot of cells for counting. Appropriately dilute in Trypan Blue (to assess viability) or 3% Acetic Acid with Methylene Blue (to assess nucleated cells). For most Leukapheresis Pack samples, a dilution of 1 in 100 is sufficient. Adjust the dilution if there are more than 100 cells per square of the hemocytometer. See Tips section for more details on performing cell counts with a hemocytometer.

NOTE: SepMate™ tubes are not intended for use with leukapheresis samples.

## Tips

For a protocol on performing total nucleated cell counts using a hemocytometer, refer to <https://www.stemcell.com/techtips>.

For a protocol on performing ammonium chloride lysis, refer to the Product Information Sheet for Ammonium Chloride Solution (Catalog #07800).

## Accessory Products

PRODUCT NAME	CATALOG #
Trypan Blue	07050
3% Acetic Acid with Methylene Blue	07060
Ammonium Chloride Solution	07800
Lymphoprep™	07801

PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED.

Copyright © 2020 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, Scientists Helping Scientists, and SepMate are trademarks of STEMCELL Technologies Canada Inc. Lymphoprep is a trademark of Alere Technologies. Spectra Optia is a registered trademark of Terumo BCT, Inc. All other trademarks are the property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.