

# TeSR™2

**Xeno-free, defined, feeder-free medium for maintenance of undifferentiated human ES and iPS cells**

Catalog #05860      500 mL Kit  
Catalog #05880      10 x 500 mL Kit



Scientists Helping Scientists™ | [WWW.STEMCELL.COM](http://WWW.STEMCELL.COM)

TOLL FREE PHONE 1 800 667 0322 • PHONE +1 604 877 0713

[INFO@STEMCELL.COM](mailto:INFO@STEMCELL.COM) • [TECHSUPPORT@STEMCELL.COM](mailto:TECHSUPPORT@STEMCELL.COM)

FOR GLOBAL CONTACT DETAILS VISIT OUR WEBSITE

## Product Description

TeSR™2 is a complete, xeno-free, serum-free, defined formulation designed for the feeder-free maintenance and expansion of human embryonic stem (ES) and induced pluripotent stem (iPS) cells in the undifferentiated state.<sup>1</sup> Complete TeSR™2 medium (Basal Medium + 5X Supplement + 250X Supplement) contains recombinant human basic fibroblast growth factor (rh bFGF) and recombinant human transforming growth factor  $\beta$  (rh TGF $\beta$ ). Addition of further growth factors is not required.

TeSR™2 may be used with Corning® Matrigel® hESC-Qualified Matrix (Corning Catalog #354277) or Vitronectin XF™ (Catalog #07180, a matrix developed and manufactured by Nucleus Biologics) as the culture matrix.

Each lot of TeSR™2 5X Supplement and 250X Supplement is used to prepare complete TeSR™2 medium and then performance-tested in a culture assay using human pluripotent stem cells (hPSCs).

## Product Information

PRODUCT NAME	CATALOG #	SIZE	COMPONENTS
TeSR™2	05860	500 mL	<ul style="list-style-type: none"><li>TeSR™2 Basal Medium (400 mL)</li><li>TeSR™2 5X Supplement (100 mL)</li><li>TeSR™2 250X Supplement (2 mL)</li></ul>
TeSR™2	05880	10 x 500 mL	<ul style="list-style-type: none"><li>TeSR™2 Basal Medium (10 x 400 mL)</li><li>TeSR™2 5X Supplement (10 x 100 mL)</li><li>TeSR™2 250X Supplement (10 x 2 mL)</li></ul>

## Component Storage and Stability

The following components are sold as complete kits (Catalog #05860 and #05880) and are not available for individual sale.

COMPONENT NAME	COMPONENT #	STORAGE	SHELF LIFE
TeSR™2 Basal Medium	05861	Store at 2 - 8°C.	Stable until expiry date (EXP) on label.
TeSR™2 5X Supplement	05862	Store at -20°C.	Stable until expiry date (EXP) on label.
TeSR™2 250X Supplement	05863	Store at -20°C.	Stable until expiry date (EXP) on label.

Please refer to the Safety Data Sheet (SDS) for hazard information.

## Preparation of Complete TeSR™2 Medium

Use sterile technique to prepare complete TeSR™2 medium (Basal Medium + 5X Supplement + 250X Supplement). The following example is for preparing 500 mL of complete medium. If preparing other volumes, adjust accordingly.

NOTE: Thaw supplements or complete medium at room temperature (15 - 25°C) or overnight at 2 - 8°C. Do not thaw in a 37°C water bath.

1. Thaw TeSR™2 5X Supplement and TeSR™2 250X Supplement. Mix each supplement thoroughly.

NOTE: Once thawed, use supplements immediately or aliquot and store at -20°C for up to 3 months. Do not exceed the shelf life of the supplements. After thawing the aliquoted supplements, use immediately. Do not re-freeze.

2. Add 100 mL of TeSR™2 5X Supplement and 2 mL of TeSR™2 250X Supplement to 400 mL of TeSR™2 Basal Medium. Mix thoroughly.

NOTE: If not used immediately, store complete TeSR™2 medium at 2 - 8°C for up to 2 weeks. Alternatively, aliquot and store at -20°C for up to 6 months. Do not exceed the shelf life of the individual components. After thawing the aliquoted complete medium, use immediately or store at 2 - 8°C for up to 2 weeks. Do not re-freeze.

If prepared aseptically, complete TeSR™2 medium is ready for use. If desired, the medium can be filtered using a 0.2 - 0.22 µm low protein binding polyethersulfone (PES) filter unit (e.g. Fisher 09-741-04 [0.2 µm, 250 mL]; Fisher SCGP00525 [0.22 µm, 50 mL]).

## Directions for Use

For complete instructions on maintaining human ES and iPS cells in TeSR™2, refer to the Technical Manual: Maintenance of Human Pluripotent Stem Cells in TeSR™2 (Document #10000005573), available at [www.stemcell.com](http://www.stemcell.com) or contact us to request a copy.

## Assessment of hPSCs

The following antibodies can be used to characterize hPSCs by flow cytometry or immunocytochemistry:

- Anti-Human SSEA-4 Antibody, Clone MC-813-70 (Catalog #60062)
- Anti-Human TRA-1-60 Antibody, Clone TRA-1-60R (Catalog #60064)
- Anti-Human OCT4 (OCT3) Antibody, Clone 3A2A20 (Catalog #60093)

For complete flow cytometry protocols and antibodies that can be used, refer to the Technical Manual: Maintenance of Human Pluripotent Stem Cells in mTeSR™1 (Document #10000005505), available at [www.stemcell.com](http://www.stemcell.com) or contact us to request a copy.

## Related Products

For related products, including specialized cell culture and storage media, matrices, antibodies, cytokines, and small molecules, visit [www.stemcell.com/hPSCworkflow](http://www.stemcell.com/hPSCworkflow) or contact us at [techsupport@stemcell.com](mailto:techsupport@stemcell.com).

## References

1. Ludwig TE et al. (2006) Derivation of human embryonic stem cells in defined conditions. *Nat Biotechnol* 24(2): 185–7.



This product was developed under license to intellectual property owned by WiCell™ Research Institute.

This product is sold for research use only (whether the buyer is an academic or for-profit entity) under a non-transferable, limited-use license. Purchase of this product does not include the right to sell, use or otherwise transfer this product for commercial purposes (i.e., any activity undertaken for consideration, such as use of this product for manufacturing, or resale of this product or any materials made using this product, or use of this product or any materials made using this product to provide services) or clinical use (i.e., administration of this product or any material using this product to humans) or the right to implant any material made using this product into an animal by, or in collaboration with, a for-profit entity, for purposes other than basic pre-clinical research applications (including without limitation teratoma assays) to validate the function of the cells. Purchasers who do not agree to the terms and conditions set forth above should return the product in acceptable conditions to the seller for a refund.

PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED. FOR ADDITIONAL INFORMATION ON QUALITY AT STEMCELL, REFER TO [WWW.STEMCELL.COM/COMPLIANCE](http://WWW.STEMCELL.COM/COMPLIANCE).

Copyright © 2021 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, and Scientists Helping Scientists are trademarks of STEMCELL Technologies Canada Inc. mTeSR and TeSR are trademarks of WARF. Corning and Matrigel are registered trademarks of Corning Incorporated. Vitronectin XF is developed and manufactured by Nucleus Biologics, and Vitronectin XF is a trademark of Nucleus Biologics. All other trademarks are the property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.