

PneumaCult™-NGEx Medium

Serum- and BPE-free medium for expansion of primary human bronchial, small airway, and nasal epithelial cells

Catalog #100-1505

500 mL



Scientists Helping Scientists™ | WWW.STEMCELL.COM

TOLL FREE PHONE 1 800 667 0322 • PHONE +1 604 877 0713

INFO@STEMCELL.COM • TECHSUPPORT@STEMCELL.COM

FOR GLOBAL CONTACT DETAILS VISIT OUR WEBSITE

Product Description

PneumaCult™-NGEx Medium is an optimized serum- and BPE-free cell culture medium for the improved expansion of primary human airway epithelial cells, including bronchial epithelial cells (HBECs), small airway epithelial cells (HSAECs), and nasal epithelial cells (HNECs). PneumaCult™-NGEx Medium builds upon previous formulations of PneumaCult™ media (e.g. PneumaCult™-Ex Plus) to support at least 15 passages of expansion, generating approximately 250X more HBECs in three weeks compared to other commercially available expansion media. The resulting cells display better-conserved differentiation potential and function, defined as the ability to form a pseudostratified mucociliary epithelium at the air-liquid interface (ALI) using PneumaCult™-ALI Medium (Catalog #05001) or a cuboidal epithelium using PneumaCult™-ALI-S Medium (Catalog #05050).

Product Information

The following components are sold as a complete kit (Catalog #100-1505) and are not available for individual sale.

COMPONENT NAME	COMPONENT #	SIZE	STORAGE	SHELF LIFE
PneumaCult™-NGEx Basal Medium	100-1504	490 mL	Store at 2 - 8°C.	Stable until expiry date (EXP) on label.
PneumaCult™-NGEx 50X Supplement*	100-1503	10 mL	Store at -20°C.	Stable until expiry date (EXP) on label.

*This product contains material derived from human plasma. Donors have been tested and found negative for HIV-1 and -2, hepatitis B, and hepatitis C prior to donation. However, this product should be considered potentially infectious and treated in accordance with universal handling precautions.

Materials Required but Not Included

PRODUCT NAME	CATALOG #
Hydrocortisone Stock Solution	07925
Animal Component-Free Cell Dissociation Kit <ul style="list-style-type: none">• ACF Enzymatic Dissociation Solution• ACF Enzyme Inhibition Solution	05426
Conical tube, 15 mL	e.g. 38009
D-PBS (Without Ca++ and Mg++)	37350
Trypan Blue	07050
Tissue culture-treated T-25 cm ² flask, with vented cap	e.g. Corning 353109

Preparation of Medium

Use sterile technique to prepare complete PneumaCult™-NGEx Medium (Basal Medium + 50X Supplement + Hydrocortisone Stock Solution). The following example is for preparing 500 mL of complete medium. If preparing other volumes, adjust accordingly.

1. Thaw PneumaCult™-NGEx 50X Supplement at room temperature (15 - 25°C). Mix gently by inverting the vial; do not vortex.

NOTE: A precipitate may be observed after thawing. This will not affect performance if the supplement is gently mixed.


NOTE: Once thawed, use immediately or aliquot and store at -20°C. Do not exceed the shelf life of the supplement. After thawing the aliquoted supplement, use immediately. Do not re-freeze.

2. Add 10 mL PneumaCult™-NGEx 50X Supplement and 0.5 mL Hydrocortisone Stock Solution to 490 mL PneumaCult™-NGEx Basal Medium. Mix thoroughly.

NOTE: If not used immediately, store complete PneumaCult™-NGEx Medium at 2 - 8°C for up to 4 weeks. Do not exceed the shelf life of the individual components.

3. If desired, add antibiotics and antifungals immediately before use (e.g. 50 µg/mL gentamicin and 250 ng/mL amphotericin B).

Protocol Diagram

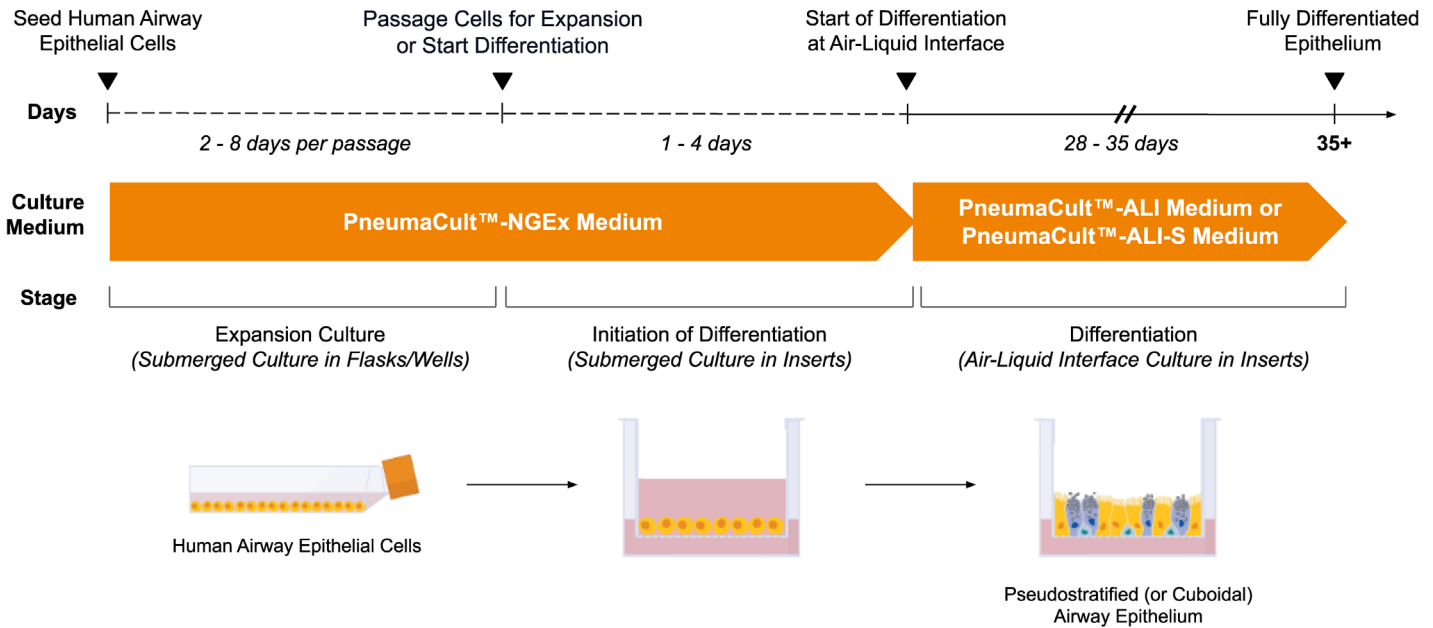

Directions for Use

Please read the entire protocol before proceeding. The following protocol is for expanding cultured (P1+) primary human airway epithelial cells (HNECs, HBECS, or HSAECs) in a single T-25 cm² flask. If using other cultureware, adjust cell numbers and volumes accordingly.

NOTE: Only use tissue culture-treated cultureware. For expanding HNECs, use tissue culture-treated cultureware coated with collagen I (e.g. Catalog #07001).

- Plate 1×10^5 cells (4×10^3 cells/cm²) in a T-25 cm² flask containing 5 mL of room temperature (15 - 25°C) complete PneumaCult™-NGEx Medium (see Preparation section).
NOTE: The optimal cell plating density may differ depending on the donor.
NOTE: If starting with cryopreserved cells, thaw cells directly into complete PneumaCult™-NGEx Medium and perform a medium change 24 hours after initial plating.
- Incubate the cells at 37°C and 5% CO₂. Perform full-medium changes every two days until cells are approximately 60 - 90% confluent and ready to be passaged. This typically takes 2 - 8 days.
NOTE: The expansion phase may take longer for some donor cell populations. To avoid weekend medium changes, perform a full-medium change with increased volume (i.e. 7.5 mL for a T-25 cm² flask) on Friday afternoon, followed by a regular full-medium change on Monday morning.
- Passage the cells using Animal Component-Free Cell Dissociation Kit, as follows:
 - Warm sufficient volumes of D-PBS (Without Ca⁺⁺ and Mg⁺⁺), complete PneumaCult™-NGEx Medium, ACF Enzymatic Dissociation Solution, and ACF Enzyme Inhibition Solution to room temperature.
 - Wash the cells once with 5 mL D-PBS. Discard the wash.
 - Add 2 mL of ACF Enzymatic Dissociation Solution and incubate the flask at 37°C for 7 - 12 minutes, or until the cells can be dislodged with gentle tapping of the flask.
NOTE: The optimal incubation time may differ depending on the donor. After 10 minutes of dissociation, continuously monitor the culture under a microscope.
NOTE: Trypsin may also be used for dissociation. However, the duration of Trypsin treatment will require optimization.
 - Add 2 mL of ACF Enzyme Inhibition Solution to the flask and transfer the cell suspension to a 15 mL conical tube.
 - Centrifuge the tube at 350 x g for 5 minutes.
 - Discard the supernatant and resuspend the cell pellet in 1 - 2 mL of complete PneumaCult™-NGEx Medium.
 - Perform a viable cell count using Trypan Blue and a hemocytometer (e.g. Catalog #100-1181).

4. To continue expansion, plate human airway epithelial cells in a new T-25 cm² flask and proceed as directed in steps 1 - 3. Alternatively, cells may be plated to initiate downstream differentiation, including air-liquid interface (ALI) cultures and airway organoids, using the following:
- PneumaCult™-ALI Medium (Catalog #05001)
 - PneumaCult™-ALI-S Medium (Catalog #05050)
 - PneumaCult™ Airway Organoid Kit (Catalog #05060)
 - PneumaCult™ Apical-Out Airway Organoid Medium (Catalog #100-0620)

For complete instructions, refer to the applicable Product Information Sheets, available at www.stemcell.com, or contact us to request a copy.

PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED. FOR ADDITIONAL INFORMATION ON QUALITY AT STEMCELL, REFER TO WWW.STEMCELL.COM/COMPLIANCE.

Copyright © 2025 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, Scientists Helping Scientists, and PneumaCult are trademarks of STEMCELL Technologies Canada Inc. All other trademarks are the property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.