



EasySep™ Human Whole Blood CD45 Depletion Kit

Positive Selection
Catalog #18289

For processing 100 mL whole blood



Scientists Helping Scientists™ | WWW.STEMCELL.COM

TOLL FREE PHONE 1 800 667 0322 • PHONE +1 604 877 0713
INFO@STEMCELL.COM • TECHSUPPORT@STEMCELL.COM
FOR GLOBAL CONTACT DETAILS VISIT OUR WEBSITE

Document #29037 | Version 1_1_1

Description

Deplete CD45+ cells directly from HetaSep™-treated whole blood.

- Fast, easy-to-use and column-free
- Up to 4 log depletion of CD45+ cells
- Isolated cells are untouched

This kit targets CD45+ cells for removal with an antibody recognizing the CD45 surface marker. Unwanted cells are labeled with antibodies and magnetic particles, and separated without columns using an EasySep™ magnet. Desired cells are simply poured off into a new tube. Isolated cells are immediately available for downstream applications such as flow cytometry, culture, or DNA/RNA extraction.

Component Descriptions

COMPONENT NAME	COMPONENT #	QUANTITY	STORAGE	SHELF LIFE	FORMAT
EasySep™ Human Whole Blood CD45 Depletion Cocktail	18289C	1 x 1 mL	Store at 2 - 8°C. Do not freeze.	Stable until expiry date (EXP) on label.	A combination of monoclonal antibodies in PBS.
EasySep™ Magnetic Nanoparticles Positive Selection	18150	1 x 1 mL	Store at 2 - 8°C. Do not freeze.	Stable until expiry date (EXP) on label.	A suspension of magnetic particles in water.
HetaSep™	07806	1 x 20 mL	Store at 15 - 25°C.	Stable until expiry date (EXP) on label.	Hetastarch solution used for erythrocyte aggregation.

PBS - phosphate-buffered saline

Components may be shipped at room temperature (15 - 25°C) but should be stored as indicated above.

Sample Preparation

For available fresh and frozen samples, see www.stemcell.com/primarycells.

WHOLE BLOOD USING HETASEP™ RED BLOOD CELL (RBC) SEDIMENTATION

1. Collect whole blood in a blood collection tube containing anticoagulant.
2. Use the minimum-sized tube for the total volume of HetaSep™ : blood sample. Add 1 part HetaSep™ to 5 parts blood and mix well.
 - For smaller samples use: 14 mL (17 x 100 mm) polystyrene round-bottom tube (e.g. Corning Catalog #352057)
 - For larger samples use: 50 mL polypropylene tubes (e.g. Corning Catalog #352070)

NOTE: For optimal results select tube sizes so that tubes can be at least 80% full, multiple tubes may be necessary. Nucleated cell recovery can decrease when using tubes less than half full due to the shorter sedimentation distance.

3. Centrifuge the sample at 50 x g for 5 minutes at room temperature (15 - 25°C) with the brake off. Remove sample from centrifuge and allow to sit undisturbed at room temperature for 5 minutes. This will allow further sedimentation of the RBCs and will improve nucleated cell recovery.

NOTE: Older samples will settle more slowly so centrifugation step at 200 x g may be required.

4. Harvest the supernatant containing the nucleated cells.
5. Top up the harvested fraction with recommended medium and centrifuge at 200 x g for 10 minutes at room temperature with the brake off.
6. Carefully remove the supernatant.
7. Resuspend cells in the recommended medium at 1/10th of the original starting volume (e.g. resuspend the cells recovered from 10 mL of whole blood in 1 mL of recommended medium).

WHOLE BLOOD USING RBC LYSIS

If RBC lysis is acceptable, lyse with Ammonium Chloride Solution (Catalog #07800).

1. Add 10 parts Ammonium Chloride Solution to 1 part whole blood.
2. Incubate for 15 minutes on ice.
3. Centrifuge at 200 x g for 10 minutes.
4. Discard supernatant, and resuspend the cells at 1/10th of the starting volume in recommended medium.

Recommended Medium

EasySep™ Buffer (Catalog #20144), RoboSep™ Buffer (Catalog #20104), or PBS containing 2% fetal bovine serum (FBS) and 1 mM EDTA. Medium should be free of Ca⁺⁺ and Mg⁺⁺.

Directions for Use – Manual EasySep™ Protocols

See page 1 for Sample Preparation and Recommended Medium. Refer to Table 1 for detailed instructions regarding the EasySep™ procedure.

Table 1. EasySep™ Human Whole Blood CD45 Depletion Kit Protocol

STEP	INSTRUCTIONS	 "The Big Easy" (Catalog #18001)
1	Prepare HetaSep™-treated sample within the volume range.	Up to 8.5 mL
	Add sample to required tube.	14 mL (17 x 100 mm) polystyrene round-bottom tube (e.g. Corning Catalog #352057)
2	Add Depletion Cocktail to sample.	100 µL/mL of sample
	Mix and incubate.	RT for 15 minutes
3	Mix Magnetic Particles. NOTE: Particles should appear evenly dispersed.	Pipette up and down more than 5 times
4	Add Magnetic Particles to sample.	100 µL/mL of sample
	Mix and incubate.	RT for 10 minutes
5	Add recommended medium to top up the sample to the indicated volume. Mix by gently pipetting up and down 2 - 3 times.	<ul style="list-style-type: none"> • Top up to 5 mL for samples < 2.5 mL • Top up to 10 mL for samples ≥ 2.5 mL
	Place the tube (without lid) into the magnet and incubate.	RT for 10 minutes
6	Pick up the magnet, and in one continuous motion invert the magnet and tube,* pouring the enriched cell suspension into a new tube.	Isolated cells are ready for use

RT - room temperature (15 - 25°C)

* Leave the magnet and tube inverted for 2 - 3 seconds, then return upright. Do not shake or blot off any drops that may remain hanging from the mouth of the tube.

Directions for Use – Fully Automated RoboSep™ Protocol

See page 1 for Sample Preparation and Recommended Medium. Refer to Table 2 for detailed instructions regarding the RoboSep™ procedure.

Table 2. RoboSep™ Human Whole Blood CD45 Depletion Kit Protocol

STEP	INSTRUCTIONS	RoboSep™ (Catalog #20000 and #21000)	
1	Prepare HetaSep™-treated sample within the volume range.	0.25 - 8.5 mL	
	Add sample to required tube.	14 mL (17 x 100 mm) polystyrene round-bottom tube (e.g. Corning Catalog #352057)	
2	Select protocol.	Human CD45 WB Depletion 18289-high recovery	
3	Mix Magnetic Particles. NOTE: Particles should appear evenly dispersed.	Pipette up and down 5 times	
4	Load the carousel.	Follow on-screen prompts	
	Start the protocol.	Press the green “Run” button	
5	Unload the carousel when the run is complete.	Isolated cells are ready for use	

Notes and Tips

OPTIMIZING RECOVERY


This no-lysis kit is designed for mid-range recovery of CD45- cells, along with 2 to 4 log depletion of CD45+ cells. If higher recovery of CD45- cells is desired, particles and cocktail may be titrated.

ASSESSING PURITY

For purity assessment of residual CD45+ cells by flow cytometry use the following fluorochrome-conjugated antibody clones:

- Anti-Human CD45 Antibody Clone HI30 (Catalog #60018) or Clone 2D1 (Catalog #60123), and
- Anti-Human CD235ab (Glycophorin A/B) Antibody, Clone HIR2 (Catalog #60111)

Data

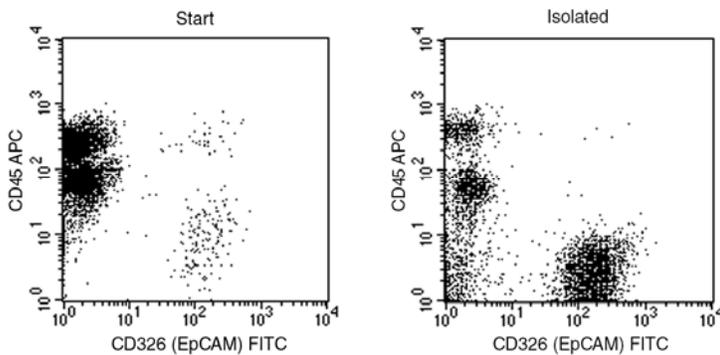

Starting with human whole blood, the depletion of CD45+ cells typically ranges from 2 - 4 log. In the example above, CAMA cells were seeded into whole blood at a starting frequency of 1.7% (98.3% CD45+). The depletion of CD45+ cells was 2 log.

NOTE: Plots were gated on Glycophorin A negative cells to exclude residual RBCs. CD326 (EpCAM) is an antibody against an epithelial cell surface antigen expressed on CAMA cells.

STEMCELL TECHNOLOGIES INC.'S QUALITY MANAGEMENT SYSTEM IS CERTIFIED TO ISO 13485. PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED.

Copyright © 2016 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, Scientists Helping Scientists, EasySep, HetaSep, and RoboSep are trademarks of STEMCELL Technologies Canada Inc. All other trademarks are the property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.