



Positive Selection  
Catalog #18167

## EasySep™ hESC-Derived CD34 Positive Selection Kit

For processing 5 x 10<sup>9</sup> cells



Scientists Helping Scientists™ | [WWW.STEMCELL.COM](http://WWW.STEMCELL.COM)

TOLL FREE PHONE 1 800 667 0322 • PHONE +1 604 877 0713  
INFO@STEMCELL.COM • TECHSUPPORT@STEMCELL.COM  
FOR GLOBAL CONTACT DETAILS VISIT OUR WEBSITE

Document #29923 | Version 1\_0\_3

## Description

Isolate highly purified CD34+ cells from human embryonic stem (ES) and induced pluripotent (iPS) cell cultures by immunomagnetic positive selection.

- Fast and easy-to-use
- Up to 99% purity
- No columns required

This kit targets CD34+ cells for positive selection with an antibody recognizing the CD34 surface marker. Desired cells are labeled with antibodies and magnetic particles, and separated without columns using an EasySep™ magnet. Unwanted cells are simply poured off, while desired cells remain in the tube. Isolated cells are immediately available for downstream applications such as flow cytometry, culture or DNA/RNA extraction.

If isolating CD34+ cells from fresh cord blood, we recommend using the EasySep™ Human Cord Blood CD34 Selection Kit II (Catalog #17896).

If isolating CD34+ cells from fresh whole blood or buffy coat, we recommend using the EasySep™ Whole Blood CD34 Selection Kit (Catalog #18086).

If isolating CD34+ cells from any other sample type, we recommend using the EasySep™ Human CD34 Positive Selection Kit (Catalog #18056).

## Component Descriptions

COMPONENT NAME	COMPONENT #	QUANTITY	STORAGE	SHELF LIFE	FORMAT
EasySep™ hESC-Derived CD34 Positive Selection Cocktail	18167C	1 x 1 mL	Store at 2 - 8°C. Do not freeze.	Stable until expiry date (EXP) on label.	A combination of monoclonal antibodies in PBS.
EasySep™ Magnetic Nanoparticles Positive Selection	18150	1 x 1 mL	Store at 2 - 8°C. Do not freeze.	Stable until expiry date (EXP) on label.	A suspension of magnetic particles in water.

PBS - phosphate-buffered saline

Components may be shipped at room temperature (15 - 25°C) but should be stored as indicated above.

## Sample Preparation

### ES or iPS CELL CULTURES

From differentiation cultures containing embryoid bodies (EBs):

1. Transfer the EBs to a 15 mL conical tube and allow to sediment.
2. Remove the medium and add 1 mL of ACCUTASE™ (Catalog #07920).
3. Incubate for 10 minutes at 37°C, then gently pipette up and down 15 times using a P1000.
4. Add 10 mL of PBS containing 2% fetal bovine serum (FBS) (Catalog #07905) and centrifuge.
5. Discard the supernatant, and resuspend the cell pellet in recommended medium.

From differentiation protocols using co-culture on adherent stromal layers:

1. Remove the non-adherent cells and transfer to a 15 mL conical tube.
2. Add 1 mL of ACCUTASE™ to the plate and incubate for 10 minutes at 37°C.
3. Remove the dissociated cells from the plate by pipetting and transfer to the same 15 mL conical tube.
4. Wash the plate with 10 mL of PBS containing 2% FBS and transfer to the same tube.
5. Centrifuge, discard the supernatant, and resuspend the pellet in recommended medium.

If clumps remain, incubate the cells with DNase I Solution (Catalog #07900) at a concentration of 100 µg/mL at room temperature (15 - 25°C) for at least 15 minutes prior to labeling and separation. Ensure the buffer used to incubate the DNase I Solution and cell suspension is free of EDTA. Filter clumpy suspensions through a 40 µm Cell Strainer (Catalog #27305) for optimal results.

After preparation, resuspend cells at the concentration indicated in Table 1 in recommended medium.

## Recommended Medium

EasySep™ Buffer (Catalog #20144), or PBS containing 2% FBS and 1 mM EDTA. Medium should be free of Ca<sup>++</sup> and Mg<sup>++</sup>.

\*\* Leave the magnet and tube inverted for 2 - 3 seconds, then return upright. Do not shake or blot off any drops that may remain hanging from the mouth of the tube.

## Notes and Tips


### ASSESSING PURITY

The EasySep™ Human CD34 Positive Selection Cocktail uses a class II anti-CD34 clone and may block some class I and II anti-CD34 clones used to assess purity by flow cytometry. Alternate clones for CD34 positive selection are available as custom kits. Contact us at [techsupport@stemcell.com](mailto:techsupport@stemcell.com) for more information. For purity assessment of CD34+ cells by flow cytometry use one of the following fluorochrome-conjugated antibody clones:

- Anti-Human CD34 Antibody, Clone 581 (Catalog #60013), Anti-Human CD34 Antibody, Clone 8G12 (Catalog #60121), AC136, or Birma K3

NOTE: Flow cytometry analysis of the positively selected cells may show slightly increased side scatter relative to the start sample.

## Data

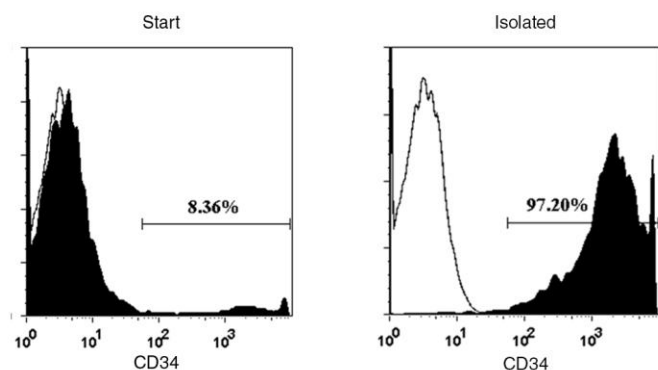

Starting with a differentiated population containing at least 5% CD34+ cells, the CD34+ cell content of the isolated fraction typically ranges from 84 - 99%. In the above example, the purities of the start and final isolated fractions are 8.36% and 97.20%, respectively.

Data courtesy of D. Kaufman, University of Minnesota.

STEMCELL TECHNOLOGIES INC.'S QUALITY MANAGEMENT SYSTEM IS CERTIFIED TO ISO 13485. PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED.

Copyright © 2016 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, Scientists Helping Scientists, and EasySep are trademarks of STEMCELL Technologies Inc. ACCUTASE is a trademark of Innovative Cell Technologies Inc. All other trademarks are the property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.