

# Dyes and Stains

## 7-AAD (7-Aminoactinomycin D)

Cell viability dye (DNA-labeling dye)



Scientists Helping Scientists™ | [WWW.STEMCELL.COM](http://WWW.STEMCELL.COM)

TOLL FREE PHONE 1 800 667 0322 • PHONE +1 604 877 0713

[INFO@STEMCELL.COM](mailto:INFO@STEMCELL.COM) • [TECHSUPPORT@STEMCELL.COM](mailto:TECHSUPPORT@STEMCELL.COM)

FOR GLOBAL CONTACT DETAILS VISIT OUR WEBSITE

Catalog # 75001.1  
75001

200 Tests  
500 Tests

## Product Description

7-AAD (7-Aminoactinomycin D) is a fluorescent cell viability dye which is excluded from live cells with intact membranes but penetrates dead or damaged cells to bind to double-stranded DNA with high affinity by intercalating between GC base pairs. It is often used as an alternative to propidium iodide (PI) to differentiate and exclude non-viable cells in flow cytometric analyses, and can be excited using 488 nm (blue), 532 nm (green), and 561 nm (yellow-green) laser lines. Compared to PI, 7-AAD offers the advantage of having minimal overlap in its emission spectra with mostly blue and green, and many red fluorophores, including the commonly used phycoerythrin (PE) and fluorescein isothiocyanate (FITC). 7-AAD is also used in DNA fluorescence imaging applications to discriminate early and late stages of apoptosis, to study cell-mediated cytotoxicity, and for chromosome banding analysis. 7-AAD has been found to exhibit antibacterial properties and exhibits growth-inhibitory activity against certain types of leukemia and sarcoma.

Chemical Name:	7-amino-actinomycin D
Alternative Names:	2,7-diamino-4,6-dimethyl-3-oxo-1-N,9-N-bis[7,11,14-trimethyl-2,5,9,12,15-pentaoxo-3,10-di(propan-2-yl)-8-oxa-1,4,11,14-tetrazabicyclo[14.3.0]nonadecan-6-yl]phenoxazine-1,9-dicarboxamide; 3H-Phenoxazine, actinomycin D deriv.; 7-Aminoactinomycin D; 7-Aminodactinomycin; Actinomycin D, 7-amino-
CAS Number:	7240-37-1
Chemical Formula:	$C_{62}H_{87}N_{13}O_{16}$
Molecular Weight:	1270.4 g/mol
Excitation Wavelength:	546 nm (DNA complex); 503 nm (free form)
Emission Wavelength:	647 nm (DNA complex); 675 nm (free form)

## Properties

Storage:	Store at 2 - 8°C. Do not freeze.
Shelf Life:	Product stable until expiry date (EXP) on label. Protect product from prolonged exposure to light.
Format/Formulation:	Phosphate-buffered saline, pH 7.2, containing 0.09% sodium azide

## Applications


Verified:	FC
Reported:	Chromosome banding, FA, FC, Fluorescence microscopy, ICC, IF
Special Applications:	This product has been verified for viability assessments of cells isolated with EasySep™ and RosetteSep™ kits.

Abbreviations: CellSep: Cell separation; ChIP: Chromatin immunoprecipitation; FA: Functional assay; FC: Flow cytometry; ICC: Immunocytochemistry; IF: Immunofluorescence microscopy; IHC: Immunohistochemistry; IP: Immunoprecipitation; RIA: Radioimmunoassay; WB: Western blotting

## Handling/Directions for Use

For flow cytometry, use 5  $\mu\text{L}$  (0.25  $\mu\text{g}$ ) of 7-AAD solution per  $1 \times 10^6$  cells in 100  $\mu\text{L}$ . Incubate cells with the dye for 5 - 10 minutes in the dark, then analyze immediately. The dye should be titrated for optimal performance for each application.

## Data/Structure

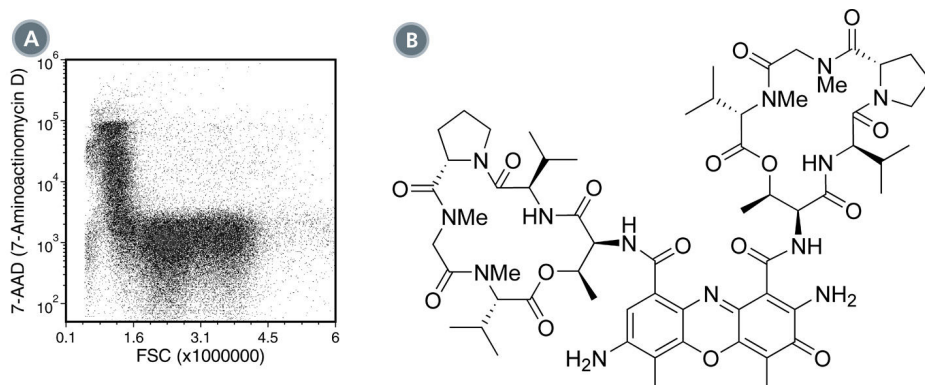

(A) Flow cytometry analysis of human peripheral blood mononuclear cells (PBMCs) labeled with 7-AAD (7-Aminoactinomycin D).  
(B) Chemical structure of 7-AAD (7-Aminoactinomycin D).

## References

1. Zembruski NCL et al. (2012) 7-Aminoactinomycin D for apoptosis staining in flowcytometry. *Anal Biochem* 429(1): 79–81. (FC)
2. Zimmermann M & Meyer N. (2011) Annexin V/7-AAD staining in keratinocytes. *Methods Mol Biol* 740: 57–63. (FC)
3. Gaforio JJ et al. (2002) Phagocytosis of apoptotic cells assessed by flow cytometry using 7-Aminoactinomycin D. *Cytometry* 49(1): 8–11. (FC)
4. Coder DM. (2001) Assessment of cell viability. *Curr Protoc Cytom* Chapter 9: Unit 9.2. (FC)
5. Lecoeur H et al. (2001) A novel flow cytometric assay for quantitation and multiparametric characterization of cell-mediated cytotoxicity. *J Immunol Methods* 253(1-2): 177–87. (FC)
6. Philpott NJ et al. (1996) The use of 7-amino actinomycin D in identifying apoptosis: simplicity of use and broad spectrum of application compared with other techniques. *Blood* 87(6): 2244–51. (FC)
7. Schmid I et al. (1994) A rapid method for measuring apoptosis and dual-color immunofluorescence by single laser flow cytometry. *J Immunol Methods* 170(2): 145–57. (FC)
8. Berg RJ et al. (1993) Flow cytometric immunofluorescence assay for quantification of cyclobutylthymine dimers in separate phases of the cell cycle. *Carcinogenesis* 14(1): 103–6. (FC)
9. Schmid I et al. (1992) Dead cell discrimination with 7-amino-actinomycin D in combination with dual color immunofluorescence in single laser flow cytometry. *Cytometry* 13(2): 204–8. (FC)
10. Toba K et al. (1992) Improved staining method for the simultaneous flow cytofluorometric analysis of DNA content, S-phase fraction, and surface phenotype using single laser instrumentation. *Cytometry* 13(1): 60–7. (FC)
11. Sengupta SK et al. (1975) 7-substituted actinomycin D analogs. Chemical and growth-inhibitory studies. *J Med Chem* 18(12): 1175–80. (FA)

## Related Products

For a complete list of related products available from STEMCELL Technologies, please visit our website at [www.stemcell.com/dyesandstains](http://www.stemcell.com/dyesandstains) or contact us at [techsupport@stemcell.com](mailto:techsupport@stemcell.com).

STEMCELL TECHNOLOGIES INC.'S QUALITY MANAGEMENT SYSTEM IS CERTIFIED TO ISO 13485. PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED.

Copyright © 2017 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, Scientists Helping Scientists, EasySep, and RosetteSep are trademarks of STEMCELL Technologies Canada Inc. All other trademarks are the property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.