

# ArciTect™ Human HPRT Positive Control Kit

**Positive control for CRISPR-Cas9 genome editing**

Catalog # 76013 1 Kit



Scientists Helping Scientists™ | [WWW.STEMCELL.COM](http://WWW.STEMCELL.COM)

TOLL FREE PHONE 1 800 667 0322 • PHONE +1 604 877 0713

[INFO@STEMCELL.COM](mailto:INFO@STEMCELL.COM) • [TECHSUPPORT@STEMCELL.COM](mailto:TECHSUPPORT@STEMCELL.COM)

FOR GLOBAL CONTACT DETAILS VISIT OUR WEBSITE

## Product Description

ArciTect™ Human HPRT Positive Control Kit is designed as a positive control for experiments using the ArciTect™ CRISPR-Cas9 genome editing system. The kit comprises ArciTect™ Human HPRT crRNA (2 nmol) and ArciTect™ Human HPRT Primer Mix (2 nmol), both of which have been tested and validated for use with the ArciTect™ line of genome editing products. HPRT, or hypoxanthine phosphoribosyltransferase, is a housekeeping gene and a commonly used control. The kit can be used to optimize transfection protocols and act as a positive control that can be used alongside custom ArciTect™ crRNAs (e.g. Catalog #76010). ArciTect™ Human HPRT crRNA first requires annealing to ArciTect™ tracrRNA (Catalog #76016) then must be combined with an ArciTect™ Cas9 Nuclease (e.g. Catalog #76001) to form a ribonucleoprotein complex. ArciTect™ Human HPRT Primer Mix can be used to amplify genomic DNA isolated from a population of transfected cells, which can subsequently be used in a T7 endonuclease assay to determine cleavage genome editing efficiency.

## Product Information

The following components are sold as a kit and are not available for individual sale.

COMPONENT NAME	COMPONENT #	SIZE	STORAGE	SHELF LIFE
ArciTect™ Human HPRT crRNA	76014	2 nmol	Store at -80°C. Alternatively, store at -20°C for up to 6 months.	Stable for 24 months from date of manufacture (MFG) on label.
ArciTect™ Human HPRT Primer Mix	76015	2 nmol	Store at -20°C.	Stable for 24 months from date of manufacture (MFG) on label.

## Materials Required But Not Included

- Nuclease-free water (e.g. Sigma Catalog # LSKNF0500)

## Directions for Use

The following instructions are for preparation of stock solutions of ArciTect™ Human HPRT crRNA (200 µM) and ArciTect™ Human HPRT Primer Mix (100 µM).

### A. ArciTect™ HUMAN HPRT crRNA 200 µM STOCK SOLUTION

1. Briefly centrifuge the vial before opening.
2. Add 10 µL of nuclease-free water. Mix thoroughly.

NOTE: If not used immediately, aliquot and store at -80°C for up to 1 month. After thawing the aliquots, use immediately. Do not re-freeze.

### B. ArciTect™ HUMAN HPRT PRIMER MIX 100 µM STOCK SOLUTION

ArciTect™ Human HPRT Primer Mix stock solution can be used in a T7 endonuclease assay to assess genome editing efficiency.

1. Briefly centrifuge the vial before opening.
2. Add 20 µL of nuclease-free water. Mix thoroughly.

NOTE: If not used immediately, aliquot and store at -80°C for up to 1 month. After thawing the aliquots, use immediately. Do not re-freeze.

For complete instructions on CRISPR-Cas9 genome editing, including annealing tracrRNA and crRNA to generate guide RNA, formation of ribonucleoprotein (RNP) complex, and transfection into target cells, refer to the Technical Bulletin: Human Pluripotent Stem Cell Genome Editing Using the ArciTect™ CRISPR-Cas9 System (Document #27084) available at [www.stemcell.com](http://www.stemcell.com) or contact us to request a copy.

## Related Products

For related products, including other genome editing tools, specialized cell culture and storage media, supplements, antibodies, cytokines, and small molecules, visit [www.stemcell.com](http://www.stemcell.com) or contact us at [techsupport@stemcell.com](mailto:techsupport@stemcell.com).

## References

- Gundry MC et al. (2016) Highly efficient genome editing of murine and human hematopoietic progenitor cells by CRISPR/Cas9. *Cell Rep* 17(5): 1453–61.
- Hultquist JF et al. (2016) A Cas9 ribonucleoprotein platform for functional genetic studies of HIV-host interactions in primary human T cells. *Cell Rep* 17(5): 1438–52.
- Kim S et al. (2014) Highly efficient RNA-guided genome editing in human cells via delivery of purified Cas9 ribonucleoproteins. *Genome Res* 24(6): 1012–9.
- Liang X et al. (2015) Rapid and highly efficient mammalian cell engineering via Cas9 protein transfection. *J Biotechnol* 208: 44–53.
- Ran FA et al. (2013) Double nicking by RNA-guided CRISPR Cas9 for enhanced genome editing specificity. *Cell* 154(6): 1380–9.
- Rupp LJ et al. (2017) CRISPR/Cas9-mediated PD-1 disruption enhances anti-tumor efficacy of human chimeric antigen receptor T cells. *Sci Rep* 7(1): 737.

STEMCELL TECHNOLOGIES INC.'S QUALITY MANAGEMENT SYSTEM IS CERTIFIED TO ISO 13485. PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED.

Copyright © 2017 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, Scientists Helping Scientists, and ArciTect are trademarks of STEMCELL Technologies Canada Inc. All other trademarks are the property of their respective holders. The purchase of the ArciTect™ products conveys to the purchaser the limited, non-transferable right to use, in accordance with applicable laws and regulations, such purchased products and any related material solely for research purposes only, not for any commercial purposes, with no warranty from the owner or assignee of the Patents, and with no liability to the owner or assignee of the Patents, or a director, trustee, officer, employee, agent, faculty, official investigator or student thereof. Distribution of ArciTect™ products is covered under at least US 8,697,359, US 8,771,945, US 8,795,965, US 8,865,406, US 8,871,445, US 8,889,356, US 8,889,418, US 8,895,308, US 8,906,616 and foreign equivalents ("Patents"). While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.