

# Human PSC-Derived Astrocytes

## Cryopreserved human PSC-derived astrocyte precursor cells for maturation into astrocytes

Catalog #70913 1 million cells XCL-1, Male



Scientists Helping Scientists™ | [WWW.STEMCELL.COM](http://WWW.STEMCELL.COM)

TOLL FREE PHONE 1 800 667 0322 • PHONE +1 604 877 0713

[INFO@STEMCELL.COM](mailto:INFO@STEMCELL.COM) • [TECHSUPPORT@STEMCELL.COM](mailto:TECHSUPPORT@STEMCELL.COM)

FOR GLOBAL CONTACT DETAILS VISIT OUR WEBSITE

## Product Description

These cryopreserved human pluripotent stem cell (PSC)-derived astrocyte precursor cells will generate a highly pure population of mature astrocytes (> 85% GFAP-positive astrocytes, < 15% class III  $\beta$ -tubulin-positive neurons) that are functional and can be maintained long-term in culture. Each vial contains  $\geq 1 \times 10^6$  viable precursor cells that can be fully differentiated and matured using STEMdiff™ Astrocyte Maturation Kit (Catalog #08550). These cryopreserved neural cells are convenient, highly consistent, and allow rapid implementation of physiologically relevant human PSC-based models for drug discovery, cell therapy validation, and neuroscientific research.

## Stability and Storage

Stable at -135°C or colder for 18 months from date of manufacture (MFG) on label.

Short-term storage of cells (< 2 weeks) at -80°C is acceptable, but should be minimized to ensure maximum stability. Upon receipt, immediately transfer vials from dry ice to storage units, avoiding exposure to room temperature. Thawed samples must be used immediately.

## Precautions

Storage of frozen cell products in the vapor phase of a liquid nitrogen storage tank is recommended. Storage in the liquid phase can result in cross-contamination if the vial breaks or is not sealed properly. Storage in the liquid phase also increases the potential for liquid nitrogen to penetrate the vial and cause it to explode when removed from storage. Use of a face shield is required as a safety precaution when transferring cells from one container to another. When handling this product do not use sharps such as needles and syringes.

STEMCELL cannot guarantee the biological function or any other properties associated with performance of cells in a researcher's individual assay or culture systems.

FOR IN VITRO RESEARCH USE ONLY. NOT APPROVED FOR DIAGNOSTIC, THERAPEUTIC, OR CLINICAL APPLICATIONS. NOT APPROVED FOR HUMAN OR VETERINARY USE IN VIVO.

## Materials Required But Not Included

| PRODUCT NAME                             | CATALOG #      |
|------------------------------------------|----------------|
| Corning® Matrigel® hESC-qualified Matrix | Corning 354277 |
| DMEM/F-12 with 15 mM HEPES               | 36254          |
| STEMdiff™ Astrocyte Maturation Kit       | 08550          |
| Falcon® Conical Tubes, 15 mL             | 38009          |
| Trypan Blue                              | 07050          |

## Directions for Use

For medium and cultureware preparation instructions (e.g. Corning® Matrigel® coating), refer to the Product Information Sheet (PIS) for STEMdiff™ Astrocyte Maturation Kit (Document #DX20345), available at [www.stemcell.com](http://www.stemcell.com) or contact us to request a copy.

### THAWING CRYOPRESERVED CELLS

NOTE: Pre-warm cultureware and media that will come in contact with cryopreserved cells. This protocol is for a single vial (1 mL) of cryopreserved cells. If using other volumes, adjust accordingly. 1 mL of cryopreserved cells can be cultured on a 60 mm cell culture dish or a 6-well plate.

1. Pre-warm two 5 mL aliquots of STEMdiff™ Astrocyte Maturation Medium at 37°C in 15 mL conical tubes.
2. Place pre-warmed tubes of medium and pre-warmed cultureware in a biosafety cabinet.
3. Remove the vial of cryopreserved cells from liquid nitrogen storage and immerse up to 2/3 of vial height in a 37°C water bath.
4. Remove vial from water bath when only a small piece of ice is visible; this will take approximately 1 minute for 1 mL of cells. Do not shake vial during thawing process.
5. Wipe the outside of the vial with 70% ethanol or isopropanol and immediately place inside the biosafety cabinet containing the pre-warmed medium.
6. Use a 1 mL pipettor to slowly transfer thawed cell suspension into one of the 15 mL conical tubes containing pre-warmed medium. Carefully transfer cell suspension dropwise into medium while swirling. Do not pipette cells up and down or generate bubbles, as cells are in a very fragile state.
7. Centrifuge diluted cell suspension at 300 x g for 5 minutes at room temperature (15 - 25°C).
8. Aspirate supernatant carefully, leaving behind a small amount of medium. Take care not to disturb the cell pellet.
9. Remove 1 mL of STEMdiff™ Astrocyte Maturation Medium from the second tube of pre-warmed medium and add to the cell pellet. **Gently resuspend** the cells by pipetting up and down slowly 4 - 6 times.
10. Perform a cell count using Trypan Blue (10 µL cell suspension + 10 µL Trypan Blue) and a hemocytometer.
11. Dilute cell suspension with an appropriate volume of the remaining STEMdiff™ Astrocyte Maturation Medium in order to achieve the required seeding density. Seed cells onto Corning® Matrigel®-coated cultureware and culture as described in section C of Directions For Use in the PIS for STEMdiff™ Astrocyte Maturation Kit (Document #DX20345), available at [www.stemcell.com](http://www.stemcell.com) or contact us to request a copy.

STEMCELL TECHNOLOGIES INC.'S QUALITY MANAGEMENT SYSTEM IS CERTIFIED TO ISO 13485. PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED.

Copyright © 2017 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, Scientists Helping Scientists, and STEMdiff are trademarks of STEMCELL Technologies Canada Inc. ACCUTASE is a trademark of Innovative Cell Technologies Inc., San Diego, CA. Corning, Falcon, and Matrigel are registered trademarks of Corning Incorporated. All other trademarks are the property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.