

# StemSpan™ T Cell Progenitor Differentiation Kit

**For expansion and differentiation of human CD34+ hematopoietic progenitor cells to the T cell lineage**

|                |        |
|----------------|--------|
| Catalog #09900 | 1 Kit  |
| Catalog #09910 | 10 mL  |
| Catalog #09920 | 0.5 mL |



Scientists Helping Scientists™ | [WWW.STEMCELL.COM](http://WWW.STEMCELL.COM)

TOLL FREE PHONE 1 800 667 0322 • PHONE +1 604 877 0713

[INFO@STEMCELL.COM](mailto:INFO@STEMCELL.COM) • [TECHSUPPORT@STEMCELL.COM](mailto:TECHSUPPORT@STEMCELL.COM)

FOR GLOBAL CONTACT DETAILS VISIT OUR WEBSITE

## Product Description

StemSpan™ T Cell Progenitor Differentiation Kit has been developed for the differentiation of CD34+ cord blood (CB) and bone marrow (BM) hematopoietic progenitor cells to T cell progenitor (pro-T) cells. StemSpan™ T Cell Progenitor Differentiation Kit typically promotes greater than 2000-fold expansion of CD5+CD7+ pro-T cells in cultures seeded with CD34+ human CB cells. After 3 weeks, on average 84% of the cultured cells express CD5 and CD7, indicative of pro-T cells, with a further subset expressing CD1a, a marker used to characterize CD5+CD7+CD1a+ pre-T cells.

StemSpan™ T Cell Progenitor Expansion Supplement (10X) contains a combination of recombinant human cytokines and other additives formulated to selectively promote the expansion and differentiation of CD34+ cells isolated from human CB and BM samples to T lineage cells. StemSpan™ T Cell Progenitor Expansion Supplement is intended for use in combination with StemSpan™ SFEM II medium, and on plates coated with StemSpan™ T Cell Differentiation Coating Material (100X).

## Product Information

The following products are components of StemSpan™ T Cell Progenitor Differentiation Kit (Catalog #09900) and are also available for individual sale.

| PRODUCT NAME   | CATALOG # | SIZE   | STORAGE         | SHELF LIFE  |
|--|-----------|--------|-----------------|---|
| StemSpan™ T Cell Progenitor Expansion Supplement (10X)   | 09910     | 10 mL  | Store at -20°C. | Stable for 12 months from date of manufacture (MFG) on label. |
| StemSpan™ T Cell Differentiation Coating Material (100X) | 09920     | 0.5 mL | Store at -20°C. | Stable for 12 months from date of manufacture (MFG) on label. |
| StemSpan™ SFEM II*                                       | 09605     | 100 mL | Store at -20°C. | Stable for 18 months from date of manufacture (MFG) on label. |

\*StemSpan™ SFEM II is also available in a 500 mL size (Catalog #09655).

## Preparation of Reagents and Materials

### StemSpan™ T Cell Progenitor Expansion Medium

Use sterile techniques to prepare StemSpan™ T Cell Progenitor Expansion Medium (StemSpan™ SFEM II + StemSpan™ T Cell Progenitor Expansion Supplement [10X]). The following example is for preparing 10 mL of medium. If preparing other volumes, adjust accordingly.

1. Thaw StemSpan™ SFEM II at room temperature (15 - 25°C) or overnight at 2 - 8°C. Mix thoroughly.  
NOTE: If not used immediately, aliquot into tubes and store at -20°C. Once aliquots are thawed, do not re-freeze.
2. Thaw StemSpan™ T Cell Progenitor Expansion Supplement (10X) at room temperature (15 - 25°C). Mix thoroughly.  
NOTE: If not used immediately, aliquot and store at 2 - 8°C for up to 1 month. Alternatively, store aliquots at -20°C. Do not exceed the shelf life of the supplement. After thawing aliquots, use immediately or store at 2 - 8°C for up to 1 month. Do not re-freeze.
3. Add 1 mL of Expansion Supplement to 9 mL of SFEM II. Mix thoroughly.  
NOTE: If not used immediately, store StemSpan™ T Cell Progenitor Expansion Medium at 2 - 8°C for up to 1 month. Do not freeze.

## StemSpan™ T Cell Differentiation Coating Material

Use sterile techniques to prepare StemSpan™ T Cell Differentiation Coating Material (Coating Material [100X] + phosphate-buffered saline [PBS; Catalog #37350]). The following example is for preparing 1 mL of StemSpan™ T Cell Differentiation Coating Material. If preparing other volumes, adjust accordingly.


1. Thaw StemSpan™ T Cell Differentiation Coating Material (100X) at room temperature (15 - 25°C). Mix thoroughly.

NOTE: If necessary, centrifuge vial in a microfuge for 30 seconds to collect liquid from cap.

NOTE: If not used immediately, aliquot and store at 2 - 8°C for up to 1 month. Alternatively, store aliquots at -20°C. Do not exceed the shelf life of the Coating Material (100X). After thawing aliquots, use immediately or store at 2 - 8°C for up to 1 month. Do not re-freeze.

2. Add 10 µL of Coating Material (100X) to 990 µL of PBS. Mix thoroughly. Use immediately.

## Protocol Diagram

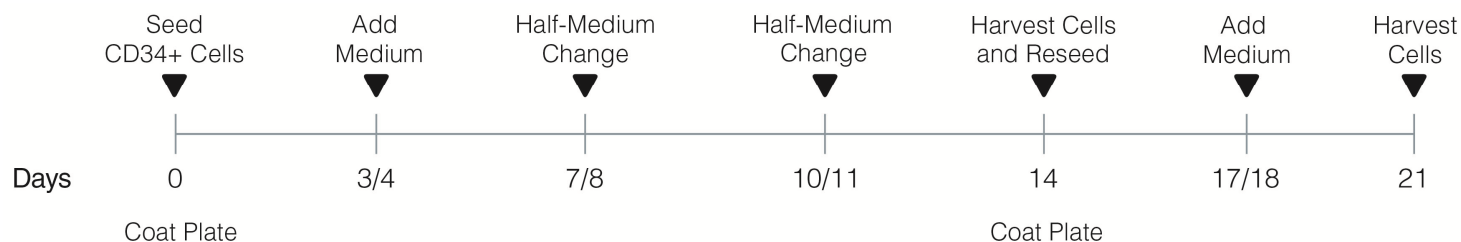

## Directions for Use

Please read the entire protocol before proceeding. Use non-tissue culture-treated cultureware.

The following instructions are for one well of a 96-well plate. If using alternative cultureware, refer to Table 1 and adjust cell numbers and volumes accordingly.

### Day 0

1. Add 100 µL of StemSpan™ T Cell Differentiation Coating Material (see Preparation of Reagents and Materials) per well of a non-tissue culture-treated 96-well plate. Refer to Table 1 for volumes required for other types of cultureware.

**Table 1. Recommended Volumes of Coating Material and Medium for Various Cultureware**

| NON-TISSUE CULTURE-TREATED CULTUREWARE | VOLUME OF STEMSPAN™ T CELL DIFFERENTIATION COATING MATERIAL | VOLUME OF STEMSPAN™ T CELL PROGENITOR EXPANSION MEDIUM | NUMBER OF CD34+ CELLS/WELL (cord blood) | NUMBER OF CD34+ CELLS/WELL (bone marrow) |
|--|---|--|---|--|
| 24-well plate                          | 500 µL/well   | 500 µL/well  | $5 \times 10^3$                         | $2.5 \times 10^4$                        |
| 12-well plate                          | 1 mL/well   | 1 mL/well  | $1 \times 10^4$                         | $5 \times 10^4$                          |
| 6-well plate                           | 2.5 mL/well   | 2.5 mL/well  | $2.5 \times 10^4$                       | $1.25 \times 10^5$                       |

2. Incubate at room temperature (15 - 25°C) for 2 hours.

NOTE: See Notes and Tips for an overnight coating method.

3. Aspirate Coating Material from the 96-well plate. Add PBS to the coated wells. Aspirate PBS just prior to use.
4. Isolate CD34+ cells from fresh (less than 72 hours old) human CB or BM using one of the following EasySep™ positive selection cell separation kits:

- Human CB : EasySep™ Human Cord Blood CD34 Positive Selection Kit II (Catalog #17896)
- Human BM: EasySep™ Human CD34 Positive Selection Kit (Catalog #18056)

NOTE: Frozen CD34+ cells from human CB or BM may also be used.

5. Perform a viable cell count using Trypan Blue (Catalog #07050) and a hemocytometer. Determine the % CD34+ cells by flow cytometry (see Notes and Tips). To determine the concentration of CD34+ cells, multiply the % CD34+ cells by the viable cell count.
6. Add purified human CD34+ cells to 100 µL of StemSpan™ T Cell Progenitor Expansion Medium (see Preparation of Reagents and Materials) as follows:
  - Human CB:  $1 \times 10^4$  CD34+ cells/mL ( $1 \times 10^3$  CD34+ cells/well)
  - Human BM:  $5 \times 10^4$  CD34+ cells/mL ( $5 \times 10^3$  CD34+ cells/well)

NOTE: This cell suspension is for one well of a 96-well plate. If using other cultureware, refer to Table 1 for volumes and cell numbers required.

7. Add 100 µL of cell suspension (prepared in step 6) to one coated well of the 96-well plate (prepared in steps 1 - 3). Incubate at 37°C.

**Day 3 or 4**

8. Carefully add 100 µL of StemSpan™ T Cell Progenitor Expansion Medium per well of the 96-well plate. Incubate at 37°C.

**Day 7 or 8**

Perform a half-medium change as follows:

9. Carefully remove 100 µL of medium from well. Do not disturb cells.
10. Add 100 µL of StemSpan™ T Cell Progenitor Expansion Medium per well. Incubate at 37°C.

**Day 10 or 11**

Perform a half-medium change as follows:

11. Carefully remove 100 µL of medium from well. Do not disturb cells.
12. Add 100 µL of StemSpan™ T Cell Progenitor Expansion Medium per well. Incubate at 37°C.

**Day 14**

13. Gently pipette cells up and down to ensure all cells are in suspension. Transfer cells to an appropriate tube.

NOTE: Harvested cells may now be used for downstream applications instead of reseeding and continuing culture for 1 week as described below. The optimal culture time is dependent on the application.

14. Coat a 96-well plate with StemSpan™ T Cell Differentiation Coating Material (see steps 1 - 3).
15. Perform a viable cell count using Trypan Blue and a hemocytometer.
16. Add cells at  $1 \times 10^5$  cells/mL to 100 µL of StemSpan™ T Cell Progenitor Expansion Medium.  
NOTE: This cell suspension is for one well of a 96-well plate. If using other cultureware, refer to Table 1 for volumes required.
17. Add 100 µL of cell suspension (prepared in step 16) to one coated well of the 96-well plate prepared in step 14 ( $1 \times 10^4$  cells/well). Incubate at 37°C.

**Day 17 or 18**

18. Carefully add 100 µL of StemSpan™ T Cell Progenitor Expansion Medium per well. Incubate at 37°C.

**Day 21**

19. Gently pipette cells up and down to ensure all cells are in suspension. Transfer cells to an appropriate tube. These pro/pre-T cells are ready for assays or analysis as required.

## Notes and Tips

- When coating cultureware with StemSpan™ T Cell Differentiation Coating Material, it may be incubated at 2 - 8°C overnight instead of at room temperature for 2 hours, if desired.
- For determining % CD34+ cells prior to plating, use one of the following fluorochrome-conjugated antibodies:
  - Anti-Human CD34 Antibody, Clone 581 (Catalog #60013)
  - Anti-Human CD34 Antibody, Clone 8G12 (Catalog #60121)
- For phenotype assessment of pro/pre-T cells by flow cytometry, use the following fluorochrome-conjugated antibodies:
  - Anti-Human CD5 Antibody, Clone UCHT2 (Catalog #60082)
  - Anti-human CD7 antibody, clone CD7-6B7
  - Anti-human CD1a antibody, clone HI149

STEMCELL TECHNOLOGIES INC.'S QUALITY MANAGEMENT SYSTEM IS CERTIFIED TO ISO 13485. PRODUCTS ARE FOR RESEARCH USE ONLY AND NOT INTENDED FOR HUMAN OR ANIMAL DIAGNOSTIC OR THERAPEUTIC USES UNLESS OTHERWISE STATED.

Copyright © 2017 by STEMCELL Technologies Inc. All rights reserved including graphics and images. STEMCELL Technologies & Design, STEMCELL Shield Design, Scientists Helping Scientists, EasySep, and StemSpan are trademarks of STEMCELL Technologies Canada Inc. All other trademarks are the property of their respective holders. While STEMCELL has made all reasonable efforts to ensure that the information provided by STEMCELL and its suppliers is correct, it makes no warranties or representations as to the accuracy or completeness of such information.